



GenePharma

Genepharma  
www.genepharma.com

## 引物探针试剂稀释说明

类型:	逆转录引物	规格:	1nmol
纯化方式:	PAGE	附送:	1.5mL DEPC 水
形态:	减压离心干燥品, 制品干燥于管底, 为干粉状		
溶液配置:	为配置成 10 $\mu$ M 溶液, 需要在每一管中加入 100 $\mu$ L DEPC 水		

类型:	PCR 引物对	规格:	1nmol
纯化方式:	PAGE	附送:	1.5mL DEPC 水
形态:	减压离心干燥品, 制品干燥于管底, 为干粉状		
溶液配置:	为配置成 10 $\mu$ M 溶液, 需要在每一管中加入 200 $\mu$ L DEPC 水		

类型:	TaqMan 探针	规格:	1nmol
纯化方式:	PAGE	附送:	1.5mL DEPC 水
标记方式	5'标记 FAM 3'标记 BHQ1		
形态:	减压离心干燥品, 制品干燥于管底, 为干粉状, 需要避光保存		
溶液配置:	为配置成 10 $\mu$ M 溶液, 需要在每一管中加入 100 $\mu$ L DEPC 水		

**说明:** 由于 DNA oligo 呈很轻的干膜附在管壁上, 打开时极易散失, 所以开管子前应先离心, 然后再慢慢打开管盖, 溶解时请加足量 DEPC 水后盖上管盖, 振荡溶解;

**贮存条件:** 减压离心干燥品可于-20 $^{\circ}$ C 保存 2-3 年, 溶液可于-20 $^{\circ}$ C 保存 6 个月-1 年, 可于 4 $^{\circ}$ C 保存 1-2 个月

---

如有疑问欢迎垂询

**电话：021-51320195(上海) 0512-86668828-8042(苏州) E-  
mail:service@genepharma.com(上海) szsupport@genepharma.com(苏州)  
http://www.genepharma.com**



GenePharma

**Genepharma**  
**www.genepharma.com**

## Primer & Probe reagent dilution mannul

<b>Kind</b>	<b>RT Primer</b>	<b>Size</b>	1nmol
<b>Purification</b>	PAGE	<b>Attachment</b>	1.5mL RNase-free H2O
<b>Format</b>	Lyophilized		
<b>Dilution</b>	Add <b>100μL</b> RNase-free H2O to <b>10μM</b> RT primer dilution		

<b>Kind</b>	<b>PCR primer set</b>	<b>Size</b>	1nmol
<b>Purification</b>	PAGE	<b>Attachment</b>	1.5mL RNase-free H2O
<b>Format</b>	Lyophilized		
<b>Dilution</b>	Add <b>200μL</b> RNase-free H2O to <b>10μM</b> primer set dilution		

<b>Kind</b>	<b>TaqMan probe</b>	<b>Size</b>	1nmol
<b>Purification</b>	PAGE	<b>Attachment</b>	1.5mL RNase-free H2O
<b>fluorescence Lable</b>	5' FAM 3'BHQ1		
<b>Format</b>	Lyophilized and keep in dark		
<b>Dilution</b>	Add <b>100μL</b> RNase-free H2O to <b>10μM</b> probe dilution		

**Storage:** Lyophilized powder could be stored at -20°C for 2-3 years, dilution at -20<sup>0</sup>C for 6 months to 1 year and at 4°C for 1-2 months

**Note:** Please open the tubes carefully and centrifuge for 5 minute at 10,000–12,000×g. Add enough RNase-free H2O to the bottom of tubes, close the tubes and vortex.

### Contact us

**Phone:** 021-51320195(Shanghai) 0512-86668828-8042(Suzhou) E-

**mail:**service@genepharma.com(Shanghai)

szsupport@genepharma.com(Shanghai)

**http://www.genepharma.com**